

Technical Data Sheet

12/07/2017

DMEM High Glucose w/ Stable Glutamine w/ Sodium Pyruvate

Product code : LM-D1110

Theoretical pH: 7.3 ± 0.3

Osmolality: 331 mOsm/kg ±10%

<u>Colour:</u> Red solution

Storage conditions: 2 to 8° C

Shelf life: 24 months

Endotoxin: <1 EU/ml

Composition: Available on request

Sterility Tests - Bacteria in aerobic and anaerobic conditions

- Fungi and yeasts

Cell Growth Test

Medium tested for the ability to support cell growth with L929 cell line

Recommended Use:

Respect storage conditions of the product

- Do not use the product after its expiry date
- Store product in an area protected from light (not necessary for saline solutions).
- Manipulate the product in aseptic conditions (e.g. : under laminar air flow)
- Wear clothes adapted to the manipulation of the product to avoid contamination (e.g. : gloves, mask, hygiene cap, overall...)

2 rue du Vieux Bourg 49340 Nuaillé Tel +33 (0)5 61 97 69 69 info@biosera.com www.biosera.com L'unité de production pour BIOSERA à Nuaillé est certifiée ISO 9001



The product is intended to be used in vitro, in laboratory only. Do not use it in therapy, human or veterinary applications.

Description

Lots of modifications of Eagle's medium have been developed since the creation of the first formulation. The most used Eagle's medium is the Dulbecco's Modified Eagle's Medium (DMEM).

It is a modification of Basal Medium Eagle (BME) that contains a concentration more important of amino acids and vitamins and also supplementary components. The original formulation contained 1000 mg/l of glucose and was used to culture embryonic mouse cells. The used of 4500 mg/l of glucose in the medium show an optimal cell growth for some cell lines.

The stable Glutamine prevents the intramolecular cyclization reaction associated with solutions of L-Glutamine. This derivative is therefore stable in solution and allows the formulation of cell culture media containing L-Glutamine that may be stored at 4°C for extended periods. The dipeptide is metabolized within the cells to yield L-Glutamine plus a second amino acid.

Uses

Supplements such as antibiotics should be added as sterile supplements to the medium. Storage conditions and shelf life of supplemented product will be affected by the nature of the supplements.

Indication of Deterioration:

Medium should be clear and free of particulate and flocculent material.

Do not use if the medium is cloudy or contains precipitate.

Other evidence of deterioration may include colour change or degradation of physical or performance characteristics.